

Updating the Driver – Window XP

To update the driver software, you need to replace the older version residing on your PC with the recent version.

The procedure for updating the Driver is described below.



Plug the connector of the USB Microscope into the PC USB port.




Click the **Start button** and choose **My Computer** from the menu.


Double-click **My Computer**, and you will see the following window.

Insert the CD-ROM with the driver file into your PC, and double-click **USB Shot Ver.1.30(E)**.

(Note) It is understood here that the driver file has been stored in the CR-ROM.

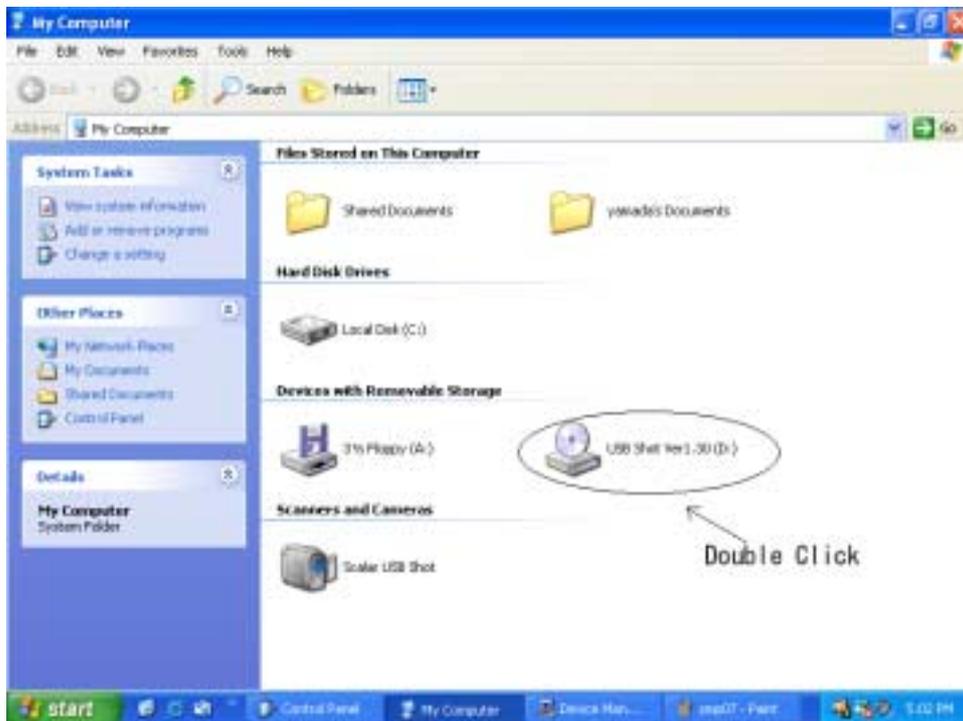

Click the folder with the **driver**

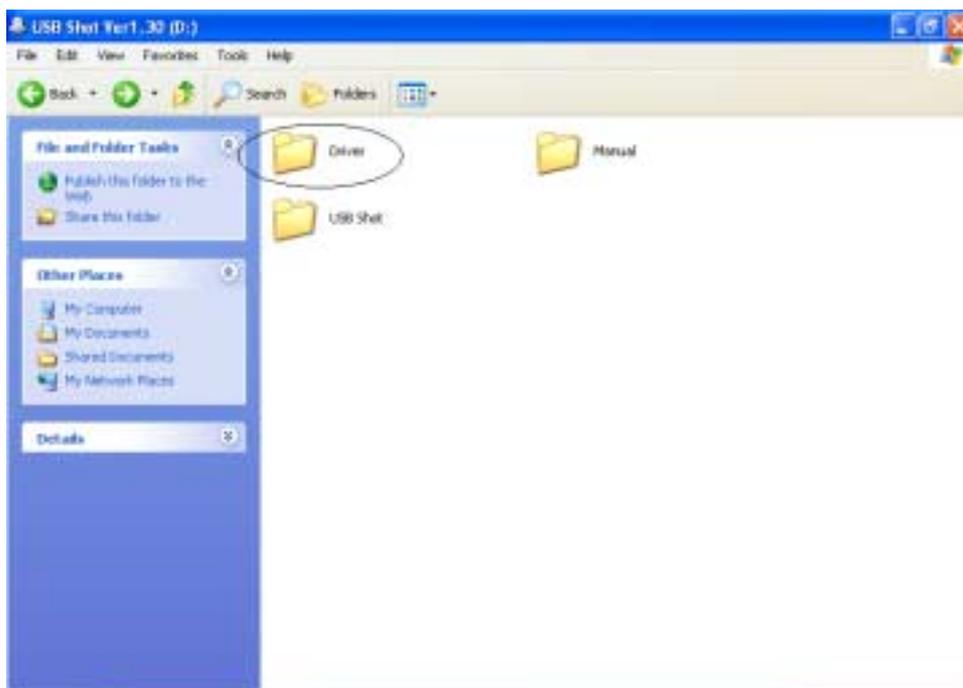




Click “Pcamrm” on the application file.



If the above message appears, click “OK” button.

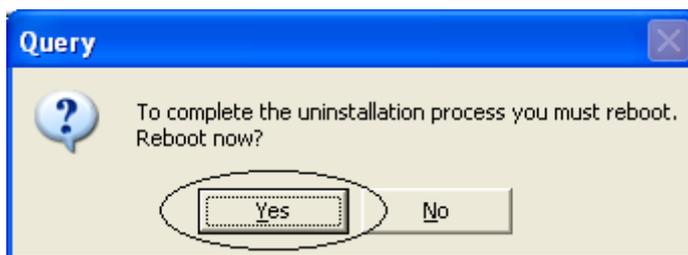

Click “Yes” button


Then rebooting begins, and so wait for some time until the PC starts up.

If you install a new driver, the following window appears when you connect the USB MicroScope with the PC. If you update the driver, the following window appears after rebooting.



Then choose installation method.

When the above window appears, it is set to “**Install the software automatically (Recommended)**” and so choose to “**Install from a list or specific location (Advanced)**” (1). Then click “**Next**”(2).

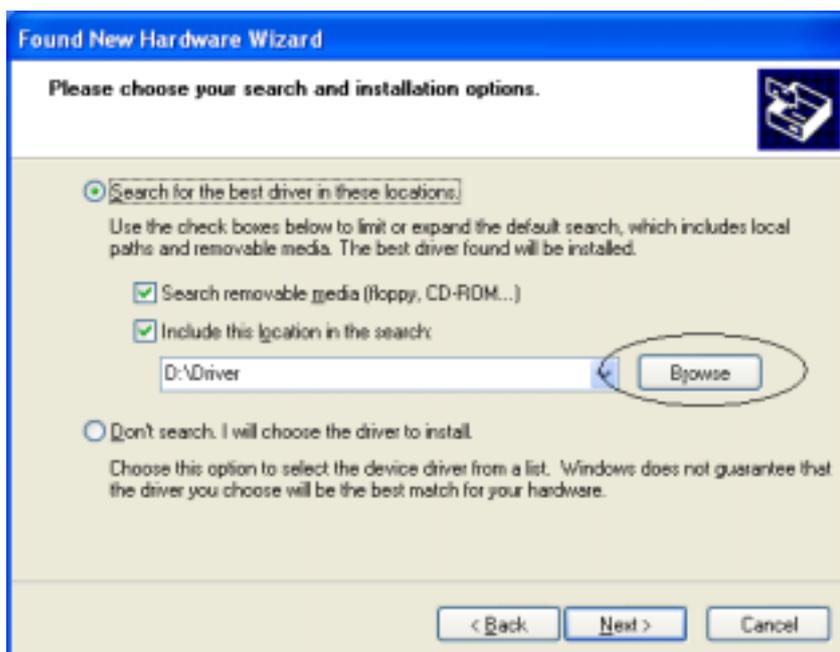


If the above window appears, click “**Browse**” button.



Designate the place where the drivers are. In this case, you choose the **driver** folder of the **CD-ROM**.

After choosing, click “**OK**” button.

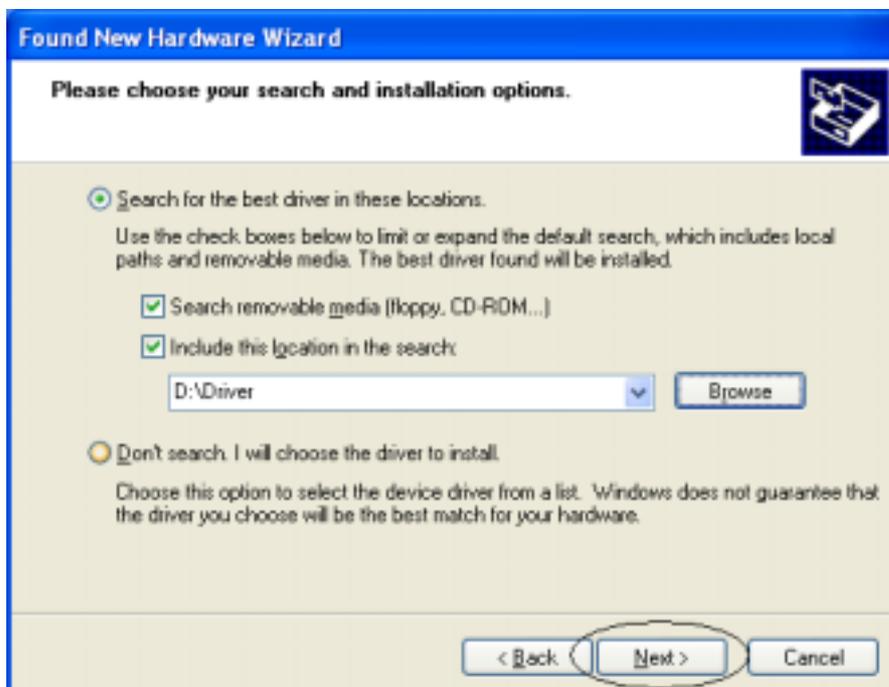



If the above window appears, click “**Next**” button.



If the above message appears, click “**Continue Anyway**” button and the installation of the driver will start.

The screen at the time of actual installation

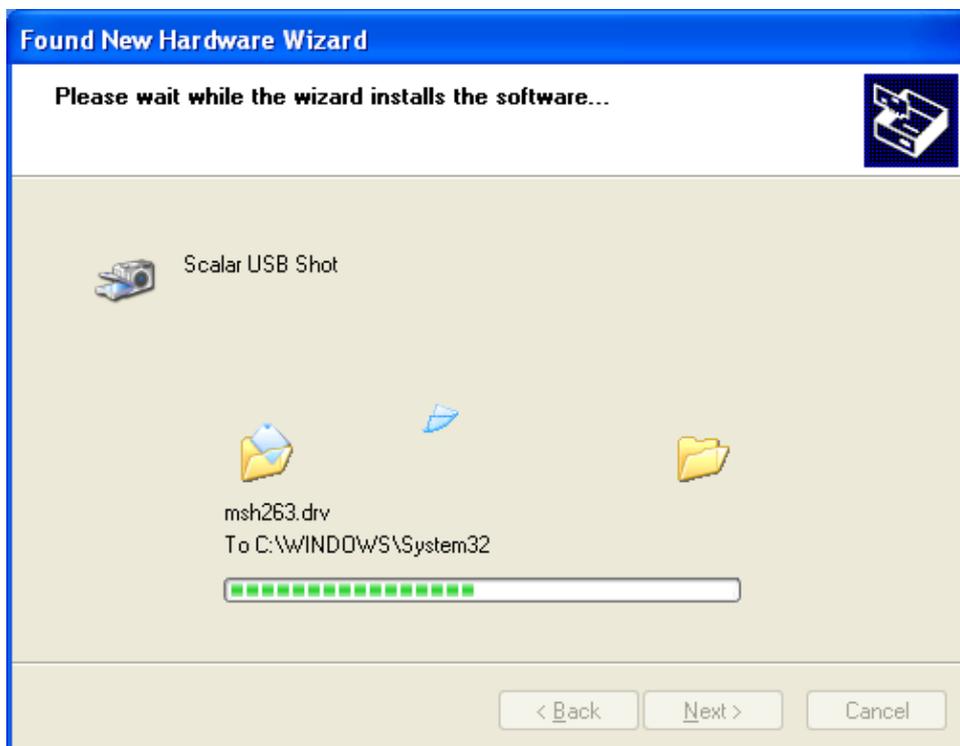

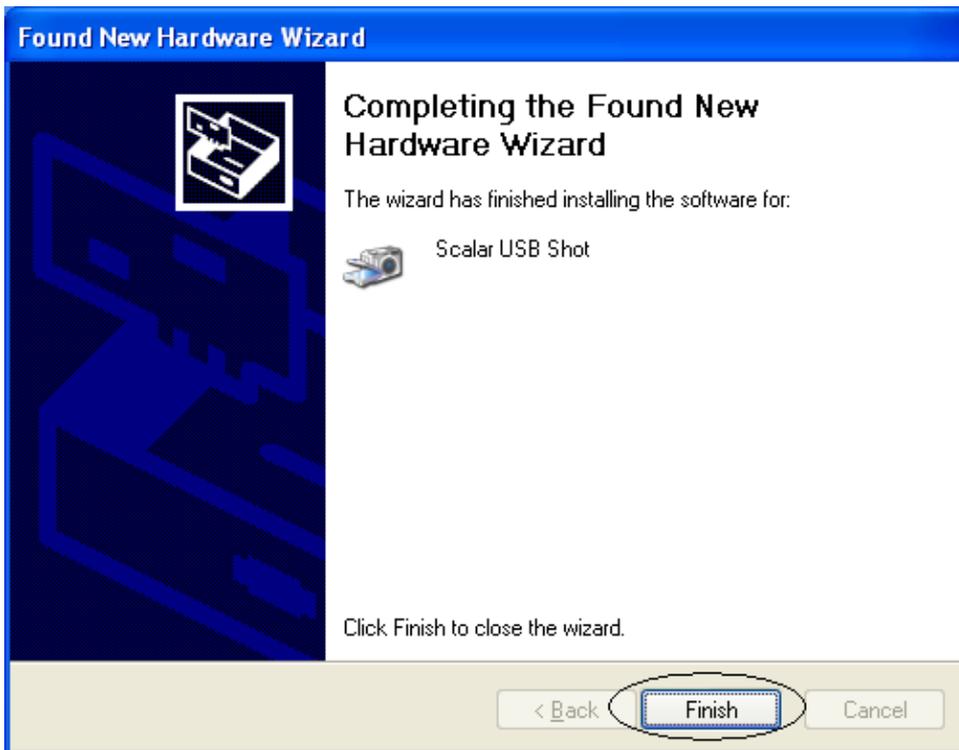

If the installation has completed, the above window will appear. Click "**Finish**" button. This completes the installation of the driver.